

A Recombinant PNGase F for the Analysis of Glycoproteins

Michael M. Rosenblatt, Chris Hosfield, Laurie Engel, Gary Kobs, and Marjeta Urh
 Promega Corporation, Madison WI
 Ravi Amunugama, David Allen, Michael Ford, and Richard Jones
 MS Bioworks LLC, Ann Arbor MI



May 21, 2012



Introduction

PNGase F is an endoglycosidase first isolated from *Flavobacterium meningosepticum*. It is a 34 kDa enzyme with specificity for N-linked glycan removal. Given the difficulty of MS/MS-based sequencing of glycopeptides, upstream removal of glycan-chains by PNGase F is a powerful tool for proteomics research.

We have expressed PNGase F as a HaloTag™ fusion and purified it to homogeneity. Recombinant PNGase F is highly active and was shown to remove N-linked glycans from a wide panel of substrates including human serum glycoproteins. Furthermore, the enzyme is active under non-denaturing conditions and therefore compatible with MS workflows.

Methods

1. Purification of Recombinant PNGase F

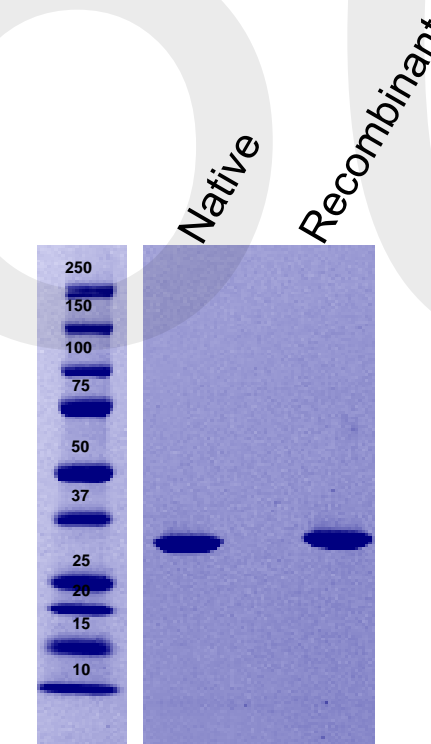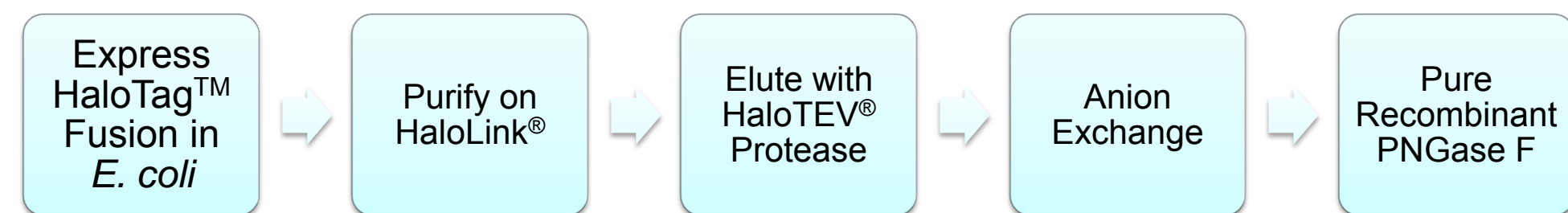


2. PNGase F Enzymatic Assays

A. Purified Glycoprotein Substrates

In a typical experiment, 50 µg of substrate was denatured by heating to 95°C for 10 minutes in 0.5% SDS + 40 mM DTT. Denatured substrate was diluted 2-fold into a final reaction buffer of 50 µL containing 1% NP-40 and 50 mM phosphate buffer, pH = 7.5. Reactions were initiated with PNGase F (500 – 2500 units) and deglycosylation was allowed to proceed for 4-18 hours at 37°C. Samples were analyzed by either SDS-PAGE (Gel-shift) or LC-MS/MS.

B. Human Serum

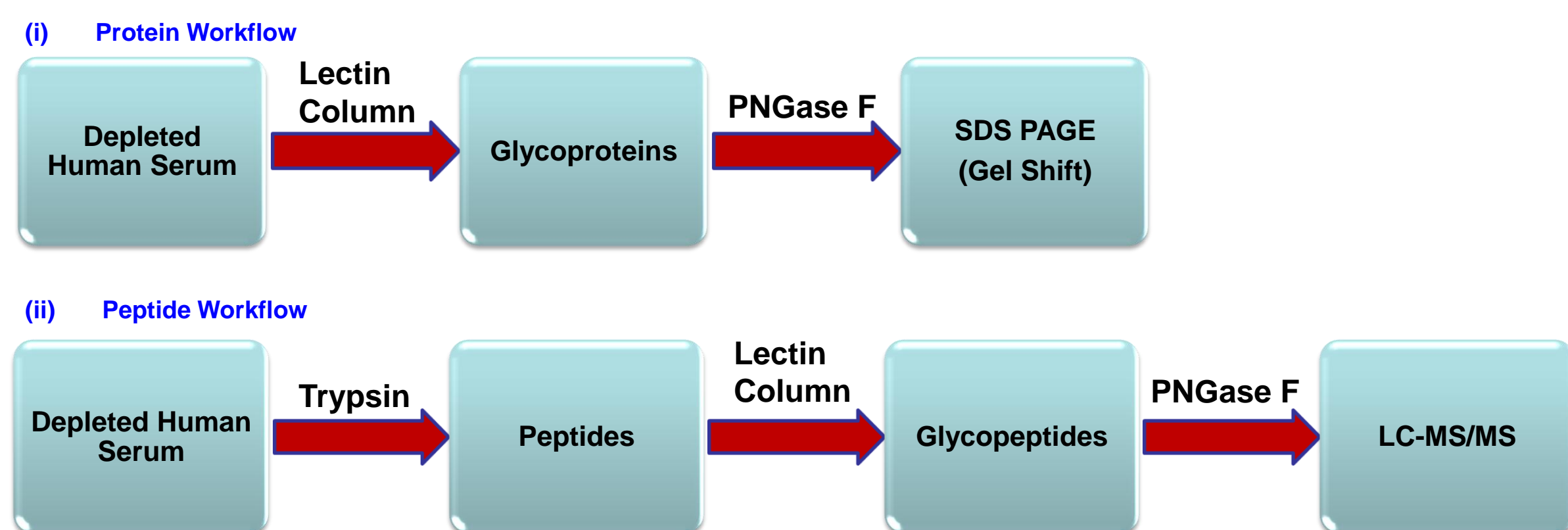

Human serum (1 mg) was depleted of albumin and IgG (Qiagen). Glycoproteins were enriched using a combined ConA/WGA lectin column. Glycopeptides were enriched by first digesting depleted human serum samples (0.2 mg) with trypsin followed by lectin enrichment. Glycopeptides were desalted with a 50 mg tC₁₈ SEP-PAK cartridge (Waters), lyophilized and reconstituted in PNGase F deglycosylation buffer. PNGase F (~1000 U) was added to the test samples while water was added to the controls.

PNGase F is Fully Active on Protein Substrates

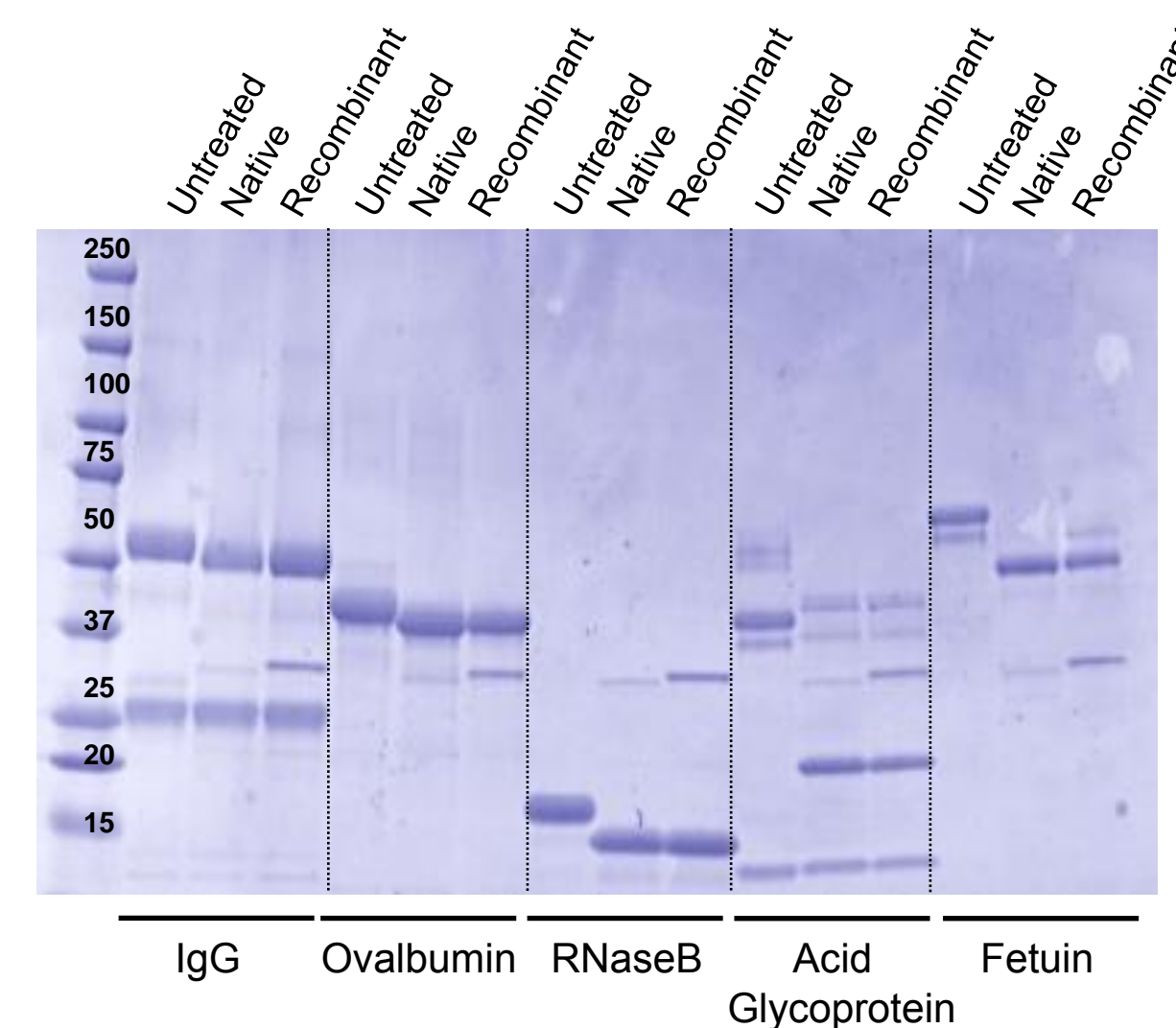


Figure 1. SDS-PAGE analysis of selected protein substrates treated with PNGase F, native and recombinant. Native and recombinant PNGase F effectively deglycosylated all substrates tested.

Immobilization of PNGase F

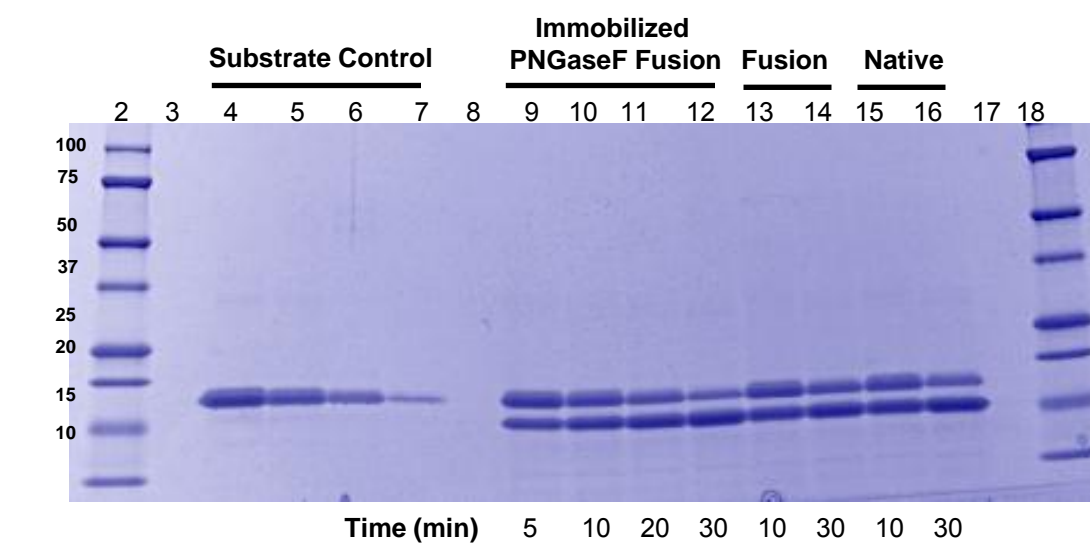


Figure 2. RNase B (100 µg) was denatured and treated with immobilized HaloTag-PNGase F, (soluble) HaloTag-PNGase F or native PNGase F at 37°C for the indicated times. Samples (5 µg) were analyzed by SDS-PAGE and visualized with SimplyBlue stain.

Analysis using LC-MS/MS

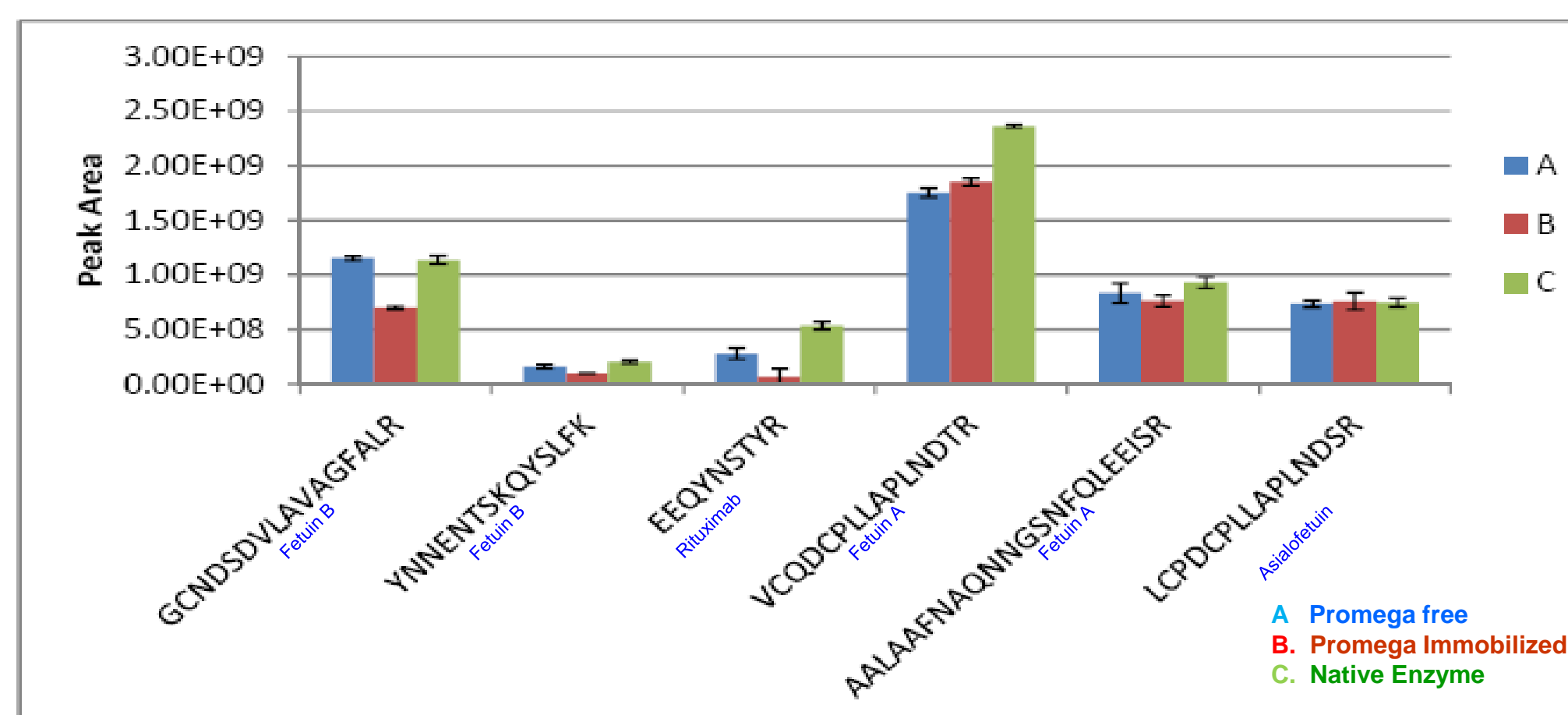


Figure 3. Specific examples of deglycosylated peptides prepared using (A) Soluble recombinant PNGase F, (B) immobilized PNGase F or (C) Native PNGase F.

PNGase F Deglycosylates Serum Proteins

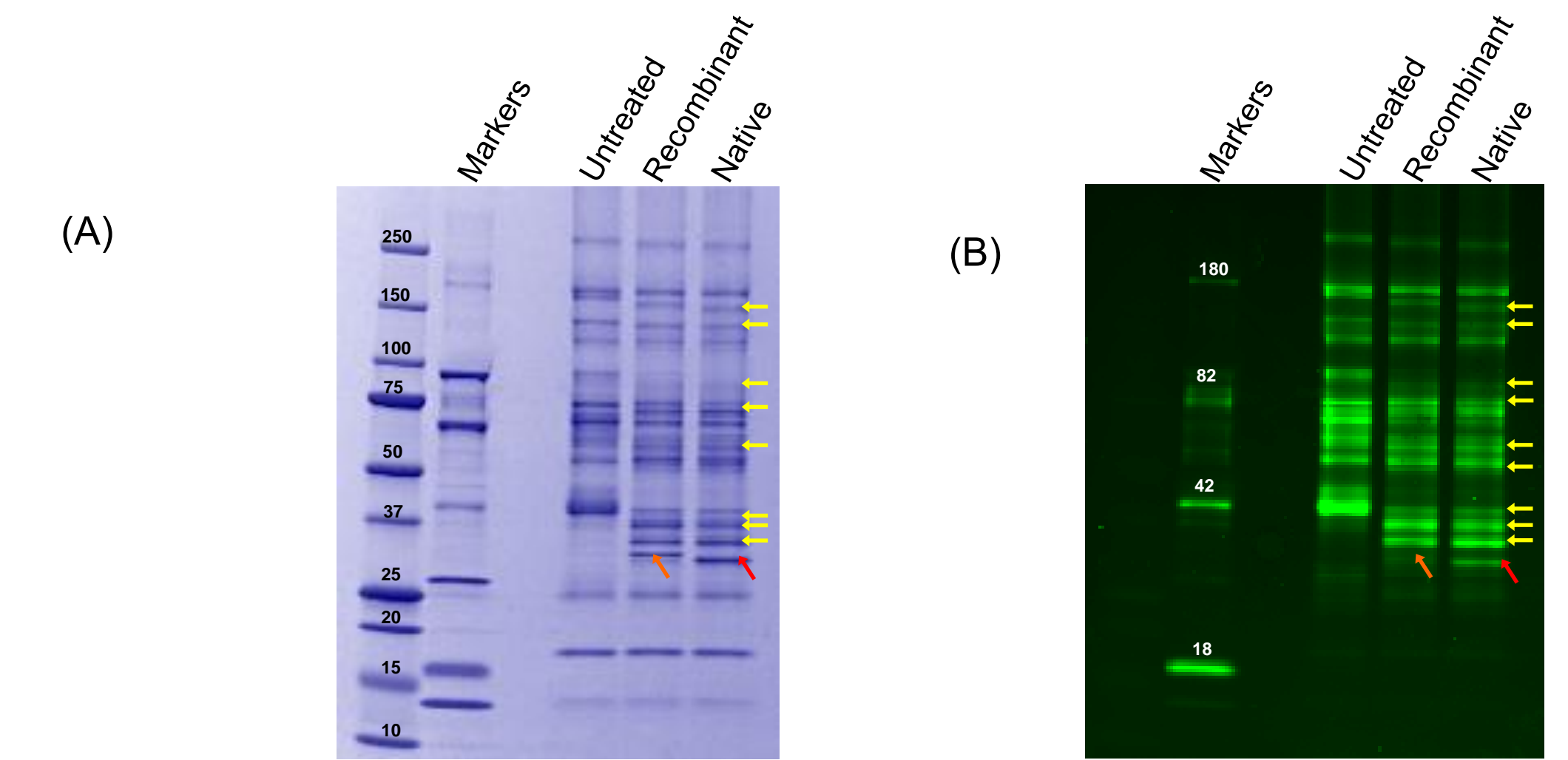


Figure 4. Enriched glycoproteins from human serum were treated with recombinant or native PNGase F at 37°C for 18 hours. Samples were analyzed by SDS-PAGE and visualized with SimplyBlue stain (A) or with Pro-Q Emerald™ glycoprotein stain (B). Arrows indicate deglycosylated proteins (yellow), recombinant (orange) and native (red) PNGase F. Deglycosylation patterns are virtually identical.

LC-MS/MS analysis of PNGase F treated Serum

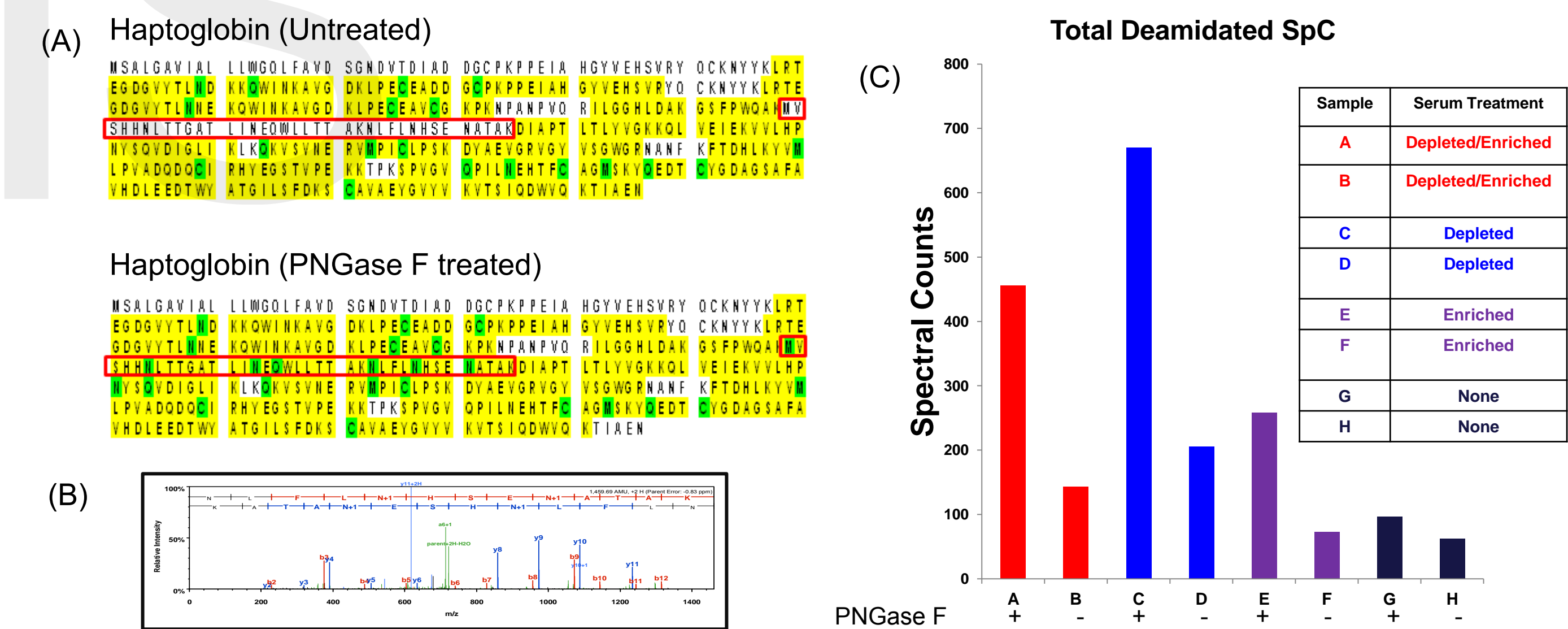


Figure 5. MS Analysis of Serum samples. (A) The sequence coverage of untreated Haptoglobin is unchanged relative to the PNGase F sample with the only exception being the identification of the glycopeptide whose spectrum is shown in panel (B). Panel (C) shows that in all cases PNGase F treatment significantly increases spectral counts for the deamidated peptides (2-5 fold)

Conclusions

- Recombinant PNGase F has been purified using HaloTag® technology to homogeneity.
- Recombinant PNGase F is fully active against both purified glycoprotein substrates and human serum glycoproteins as compared with native PNGase F.
- PNGase F improves sequence coverage is required to identify glycosylation sites.
- Deglycosylation of serum using PNGase F is compatible with MS workflows (i.e. can be used with non-denaturing buffers).
- Immobilization of PNGase F might be a useful tool for automation of sample prep workflows.